Beyond the Genome:

Unraveling Protein Variability with Quantum-Si's Next-Generation Protein SequencingTM Technology


Khanh D. Q. Nguyen, Mathivanan Chinnaraj, Sharmin Sultana, Alexa Andrzejewski, Marla Charron, Cassie Simonides, Kenneth Skinner, Michael Meyer, Abde Ali Kagalwalla, Ben Moree, and Brian Reed Quantum-Si Incorporated, 29 Business Park Drive, Branford, CT 06405

INTRODUCTION

Protein sequencing is a groundbreaking advancement in proteomics that augments genomics and transcriptomics research by providing crucial insights into the functional proteins encoded by the genome. Protein sequencing offers a more complete understanding of cellular processes and disease mechanisms by detecting changes at the protein level, such as post-translational modifications (PTMs), which cannot be captured by genomics data alone. Next-Generation Protein Sequencing[™] (NGPS) on Platinum[®] enables researchers to identify and characterize proteins with single-molecule resolution in a simple workflow and on a benchtop instrument.


Distinction of Variants of SARS-CoV-2 Virus

Kinetic signatures enable distinction of three major variants of Covid spike proteins through just two single-amino-acid substitutions.

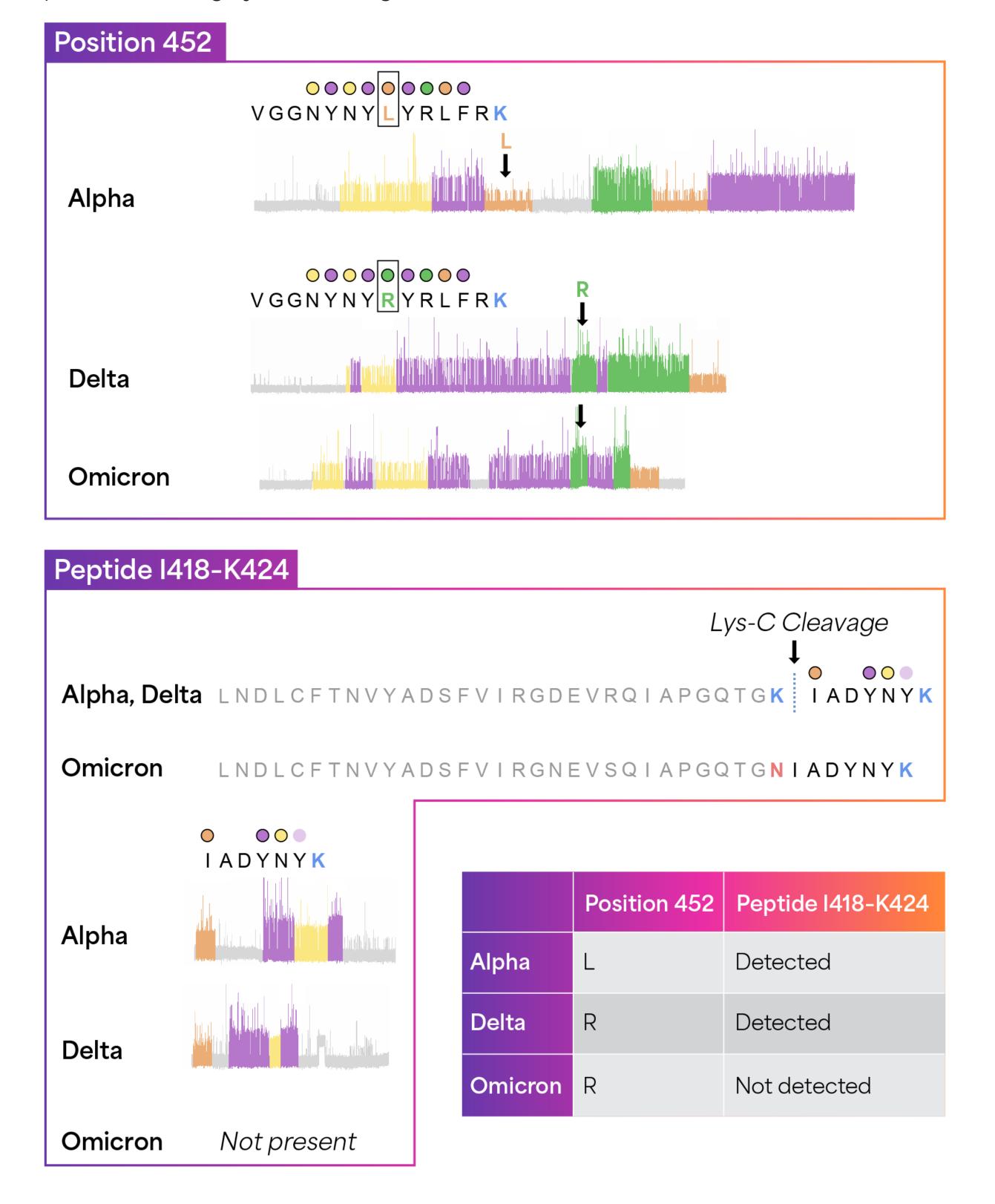

Accurate Identification of HSA from Human Serum Extracted by SDS-PAGE

The new protein inference tool (7,921-protein panel) enables accurate identification of HSA from extracted SDS-PAGE gel band of human serum.

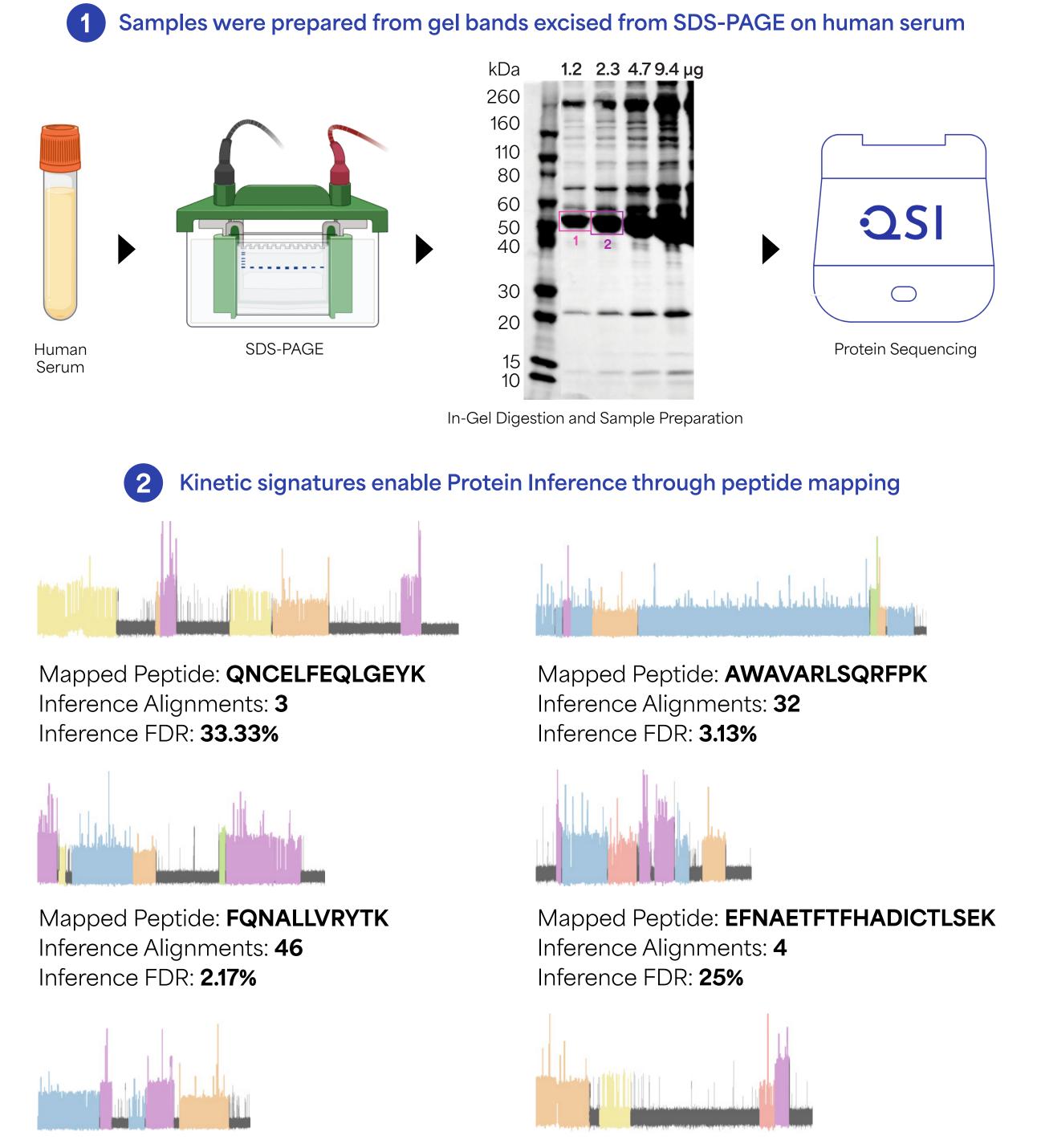

To demonstrate the versatility of Platinum and the use of kinetic signatures, we sequenced various types of samples, including protein variants with single-amino-acid changes, mixtures of recombinant proteins, peptides with PTMs, proteins immunoprecipitated from human serum, and proteins isolated from human serum via fractionation with SDS-PAGE. First, Platinum was utilized to successfully discern a single amino acid substitution at the 12th position of a peptide in ubiquitin, showcasing the sensitivity of the system in the discovery of mutations deep into peptides. Next, we **successfully dis**criminated three variants of SARS-CoV-2 spike proteins: Alpha, Delta, and **Omicron** using two single amino acid substitutions among these variants.

Next, we sequenced a mixture of ten recombinant proteins: HSA, VIME, IL6, PDL1, APOE4, FGF2, AKT1, CDNF, IL4, and H4. The resulting peptides generated distinct kinetic signatures aligned to their respective sequences, highlighting the efficacy of Quantum-Si's sequencing platform in analyzing multi-protein mixtures at reduced input concentrations. Additionally, we demonstrated the power of Platinum to detect PTMs on the basis of kinetic changes by detecting citrullination and dimethylation of arginine-two PTMs that play key roles in disease states such as cardiovascular disease, autoimmune disease, and cancer.

Finally, we developed software based on a statistical inference method to identify proteins from sequencing data without prior knowledge via mapping to a large reference panel consisting of a subset of the human proteome. This software can also map to user-specified panels, enabling protein identification tailored to specific biological pathways. To demonstrate this capability, we isolated proteins from human serum via immunoprecipitation or SDS-PAGE and correctly identified them from the sequencing data with high confidence using an 8,000-protein reference panel.

Accurate Identification of a Mixture of 10 Proteins

A mixture of 10 recombinant proteins of various molecular weights and biological functions was successfully sequenced with Platinum.

10-Protein Mix - 24,720 Alignments

Mapped Peptide: LVNEVTEFAK Inference Alignments: **10** Inference FDR: 10%

Inference Alignments: 8

Inference FDR: 12.5%


Mapped Peptide: **AVMDDFAAFVEK**


Mapped Peptide: YLYEIARRHPYFYAPELLFFAK Inference Alignments: **14** Inference FDR: 7.14%



METHODS


- Proteins are reduced, alkylated, and digested with LysC.
- Peptides are functionalized, conjugated, and immobilized on the surface of a proprietary semiconductor chip.
- Fluorescently labeled N-terminal amino acid (NAA) recognizers and aminopeptidases are added to the semiconductor chip.
- Fluorescent intensity and duration of each NAA binding event generates a unique kinetic signature.
- Kinetic signatures are analyzed to align reads to reference peptides and compute false discovery rate (FDR).

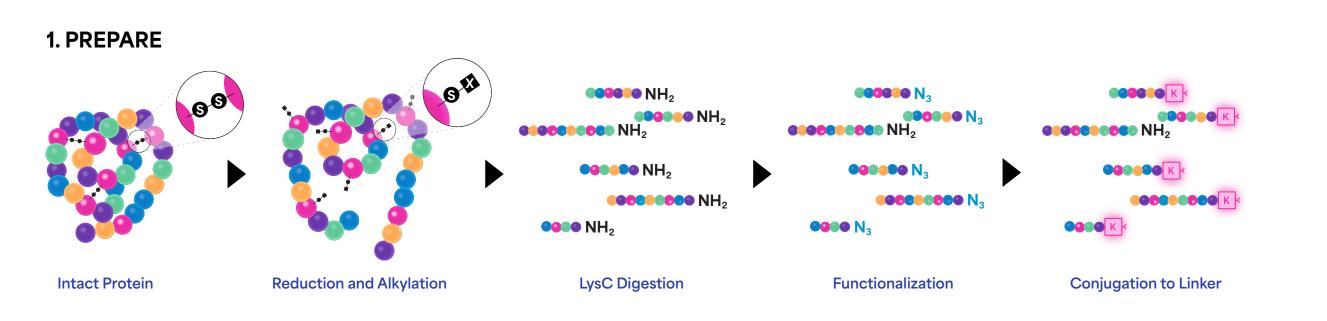

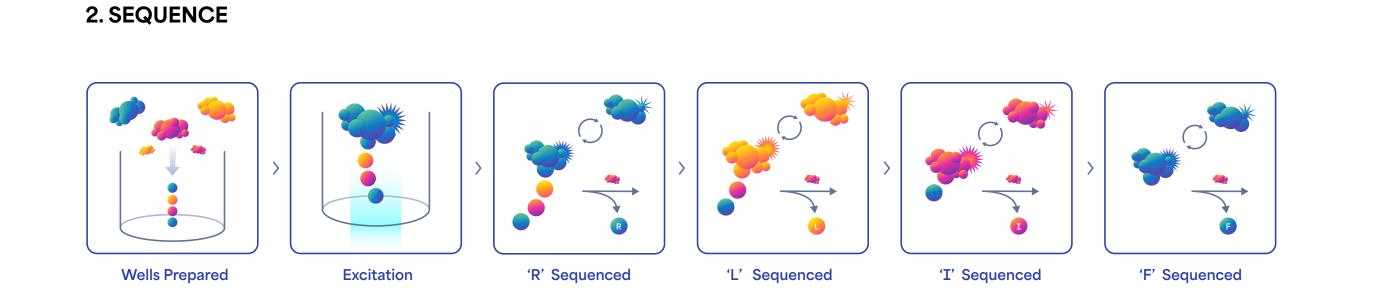

2	ANA		7 C
Э.	AINF	1LIZ	-

IL4 126 residues 22,160 Reads	EAN (103-117): 15,548 Alignments	
0 H K* 2 2 C D I T L Q E I I K* 56 12 12 T L N S L T E Q K* 27 21 T L C T E L T V T D I F A A S K* 4 37	Coverage F A N Q S T L E N F L 97.1% 38.8% 51.3% 65.0% 3.5% 97.4% 47.5% 39.5% 72.1% 77.5%	
N T T E K* 42 42 E T F C R A A T V L R Q F Y S H H E K* 61 61 D T R C L G A T A Q Q F H R H K* 77	PD (s) 0.33 0.17 0.41 1.01 0.22 0.77 0.29 0.42 2.35 1.19	
Q L I R F L K ⁴⁶⁴¹ R L D R N L W G L A G L N S C P V K ⁴⁶¹ 1405 1405 E A N Q S T L E N F L E R L K ⁴ 117	IPD (s) 11.8 15.8 11.5 7.12 17.0 4.69 18.1 14.2 4.49 6.39	
T I M R E K* 123 123 Y S K* 126	ROI Start (m) 2.86 31.8 57.8 81.1 78.4 99.7 138.0 174.0 186.0 201.0 217.0	
117 123 123 126 126 ★ Alignment Count	ROI Duration (m) 17.8 34.3 19.6 15.6 30.9 24.1 51.3 22.6 12.6 8.55 96.0	

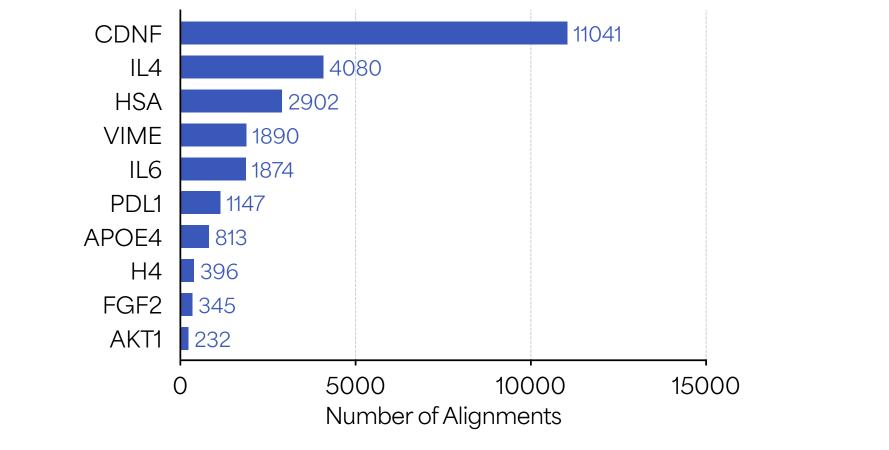


PD (s)


Example Trace

Coverage

PD (s)

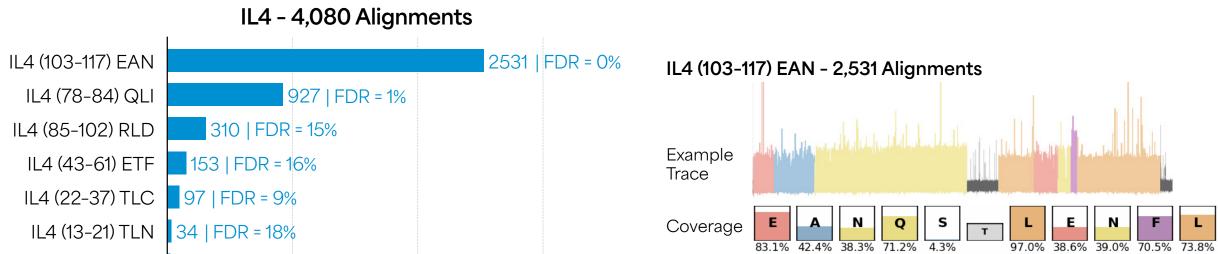


3000 1000 2000 Number of Alignments

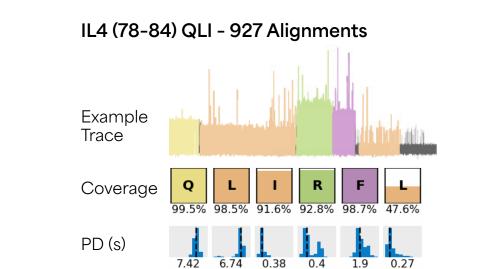

8 | FDR = 32%

IL4 (3-12) CDI 28

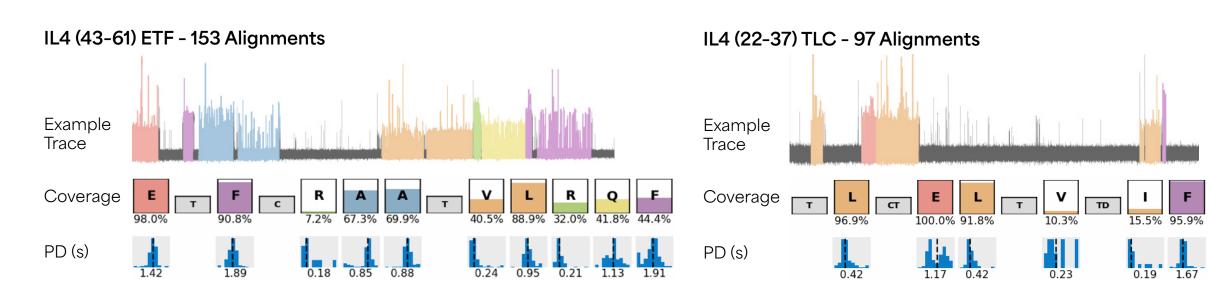

IL4 (13-21) TLN - 34 Alignments

IL4 (3-12) CDI - 28 Alignments



L P R N L W G L A 99.7% 8.4% 98.1% 97.1% 16.5% 23.9% 16.5%

IL4 (85-102) RLD - 310 Alignments


0.81 0.3 0.45 2.49 1.27

30 min

3 ALBU has a probability of 99.99% to be present in the sample


Protein Inference - 99.99% Probability for ALBU

Peptide Alignment for ALBU

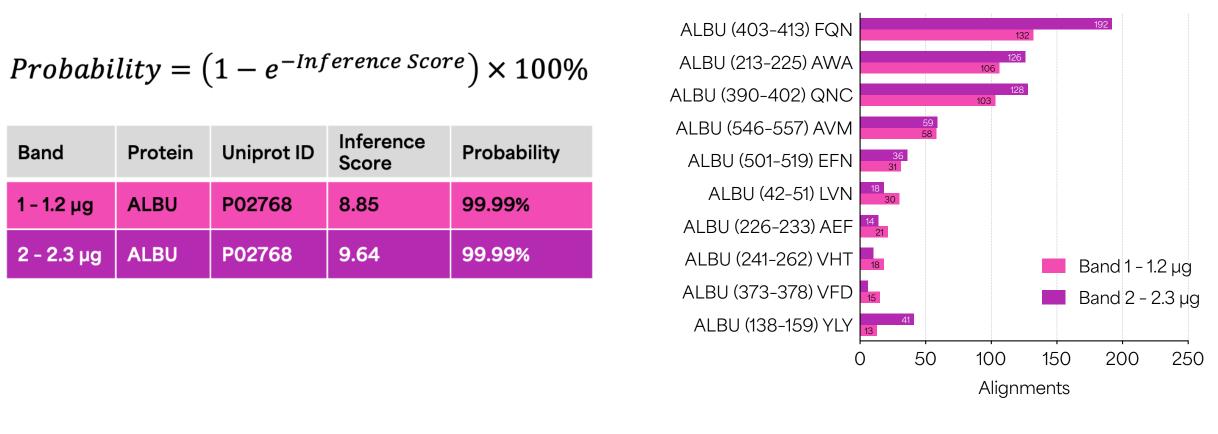


Accurate Identification of IL6 Immunoprecipitated from Human Serum


IL6 immunoprecipitated from human serum was correctly identified as the top protein against an 7,921-protein reference panel.

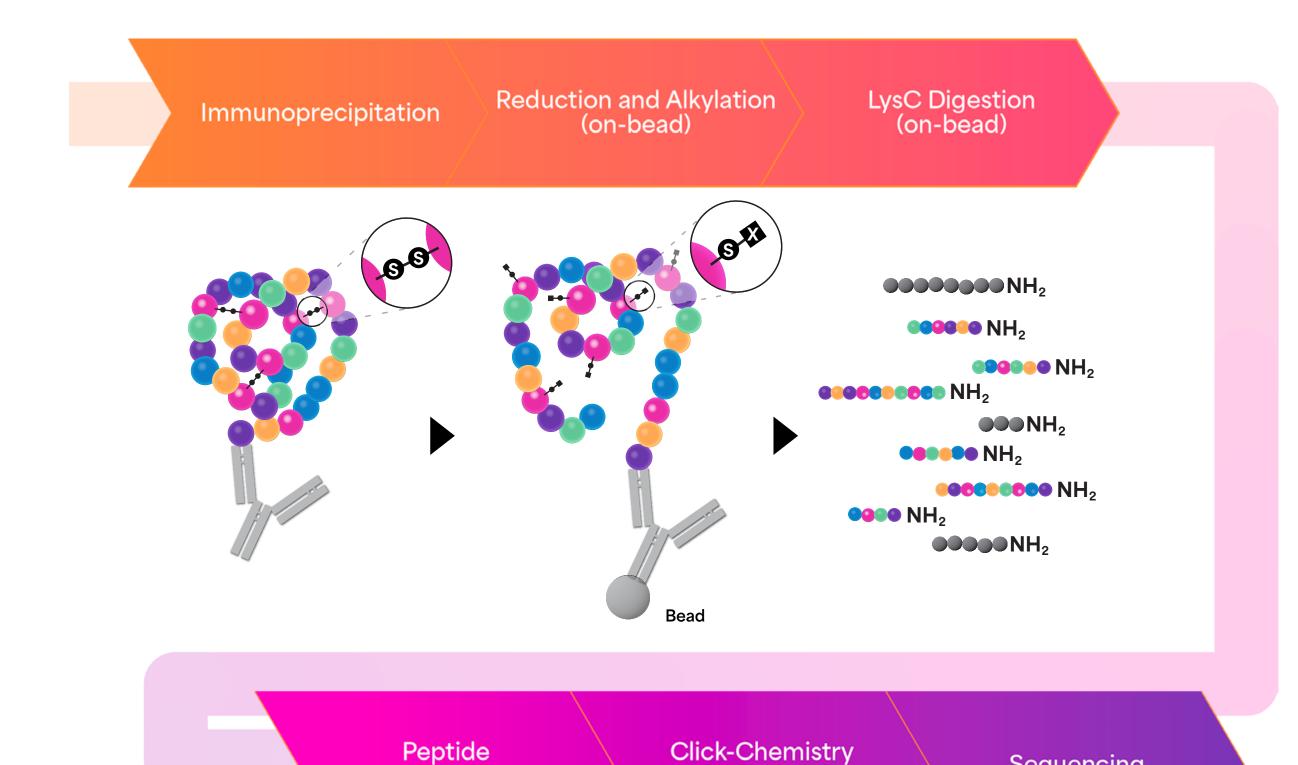

RESULTS

Detection of a Single Amino Acid Variant at the 12th Position of a Ubiquitin Peptide

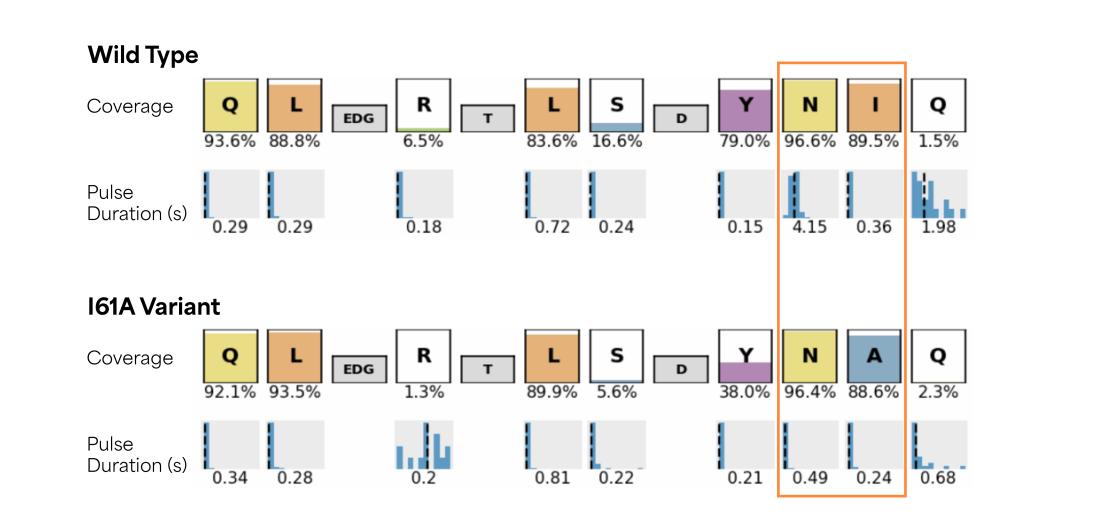


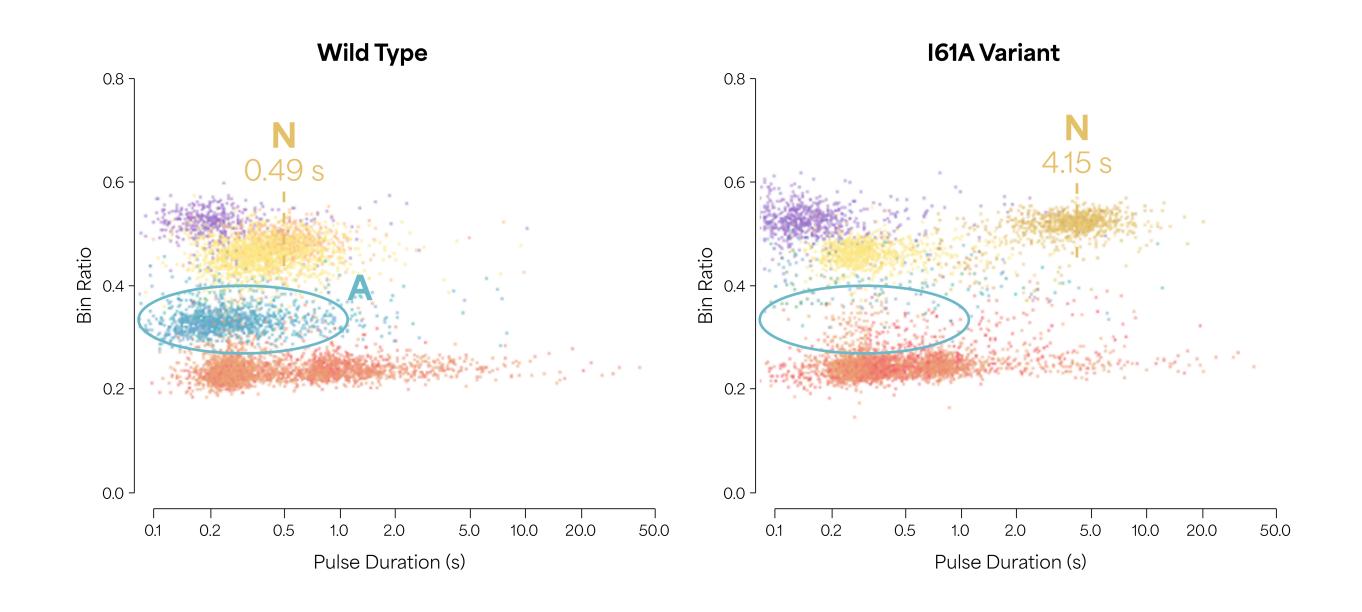




Detection of Arginine Citrullination from Peptides Derived from Vimentin

Citrullination led to the recognition of the first V residue, while the recognition of the second R residue was eliminated.

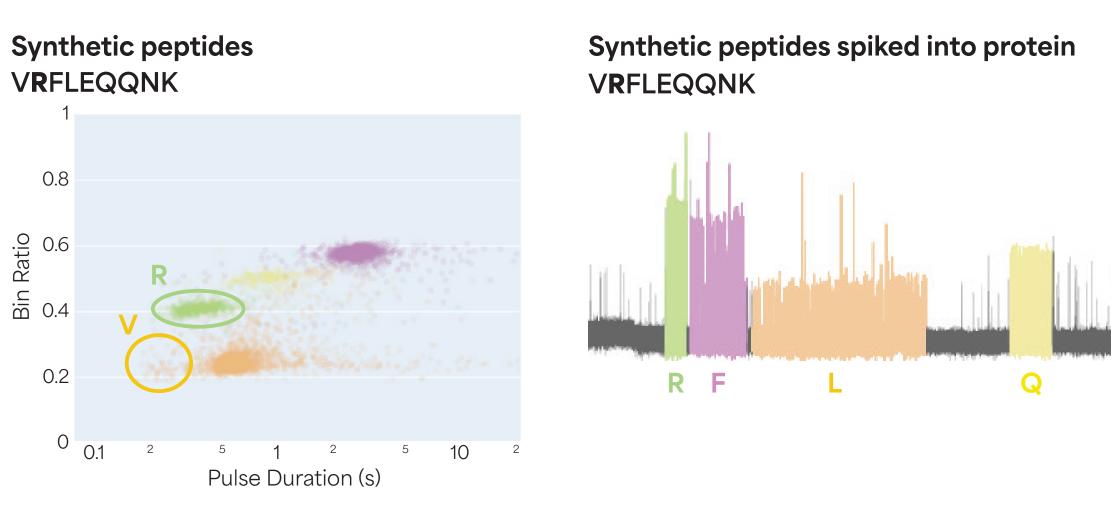


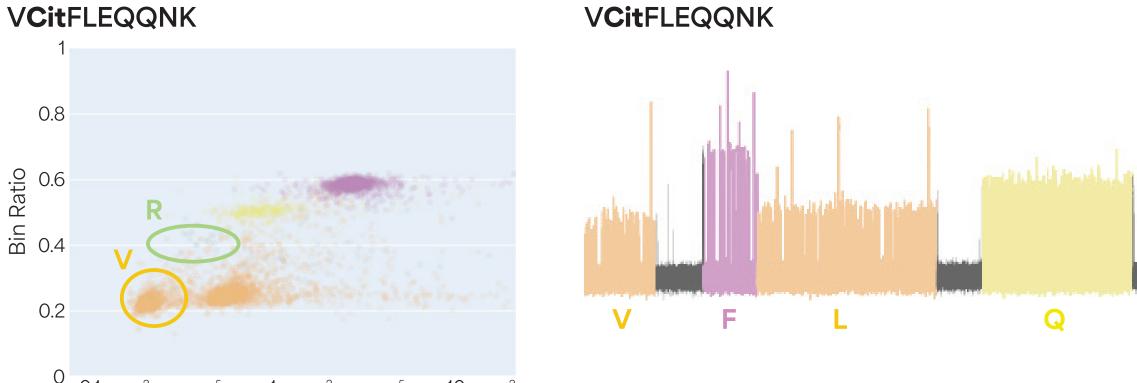

0 0.1 ² ⁵ 1 ² ⁵ 10 ² Pulse Duration (s)


0.8

0.2

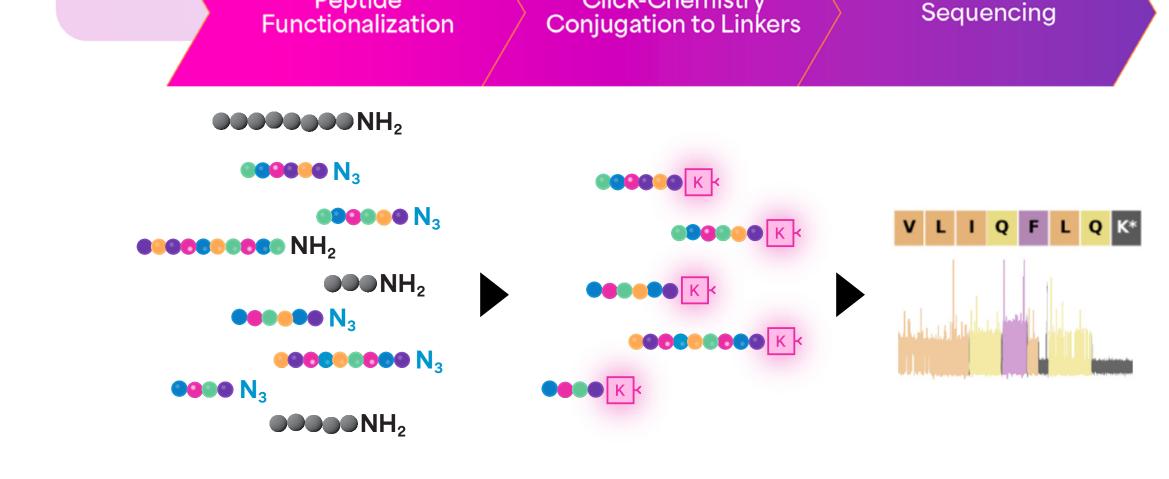

Protein Inference - 99.97% Probability for IL6

Peptide Alignment - 5,803 Alignments for IL6

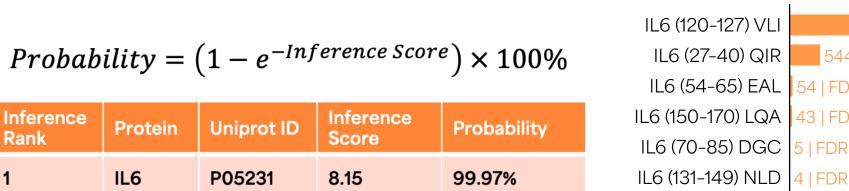

5149 | FDR = 0% IL6 (27-40) QIR 544 | FDR = 2% 54 | FDR = 4% 13 | FDR = 9% | FDR = 40% 4 | FDR = 75% IL6 (89-119) IIT I FDR = 100% 6000 Alignments

REFERENCE

Brian D. Reed et al, Science 2022, 378 (6166) 186-192.

TRADEMARKS/LICENSING

All trademarks are the property of Quantum-Si, Inc. or their respective owners. For specific trademark information, see: www.quantum-si.com/privacy-policy.

For research use only. Not for use in diagnostic procedures.